Workshop BWAW

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

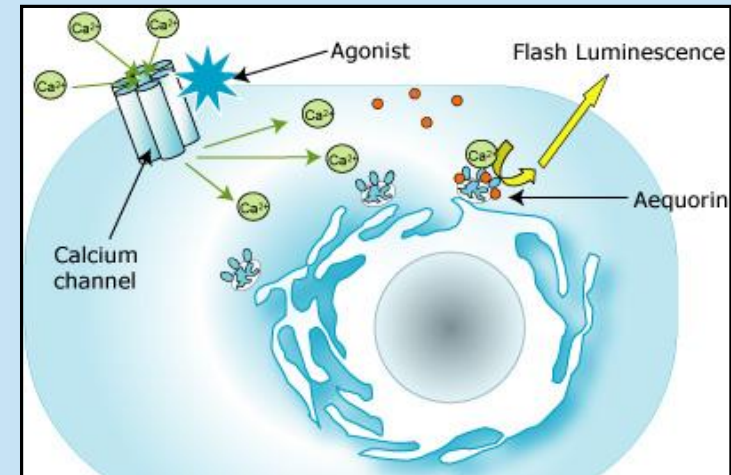
Liquid Handling in High Throughput Screening (HTS)

Highly **precise** and **parallel** pipetting and dispensing of **smallest** liquid volumes into microplates are fundamental processes in high throughput biological und chemical analysis.

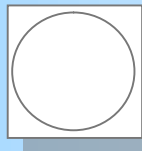
Microplate formats


96 wells

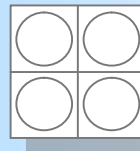
384 wells

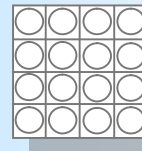1536 wells

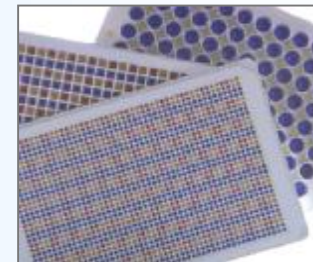10,000 biological tests need

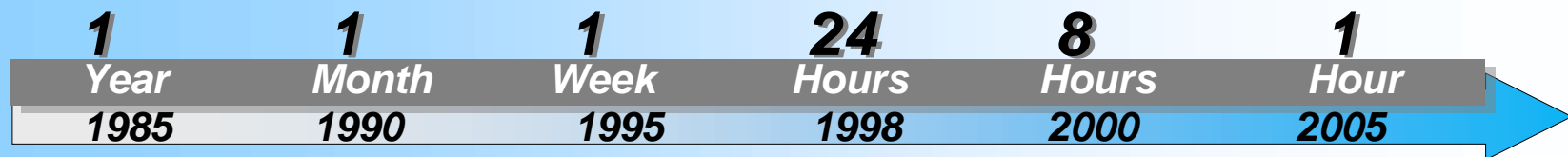Core Competence

more than 20 years experience in development

CyBi[®]-NanoJet

Sumal MD 96

“single-arm bandit”

Liquid Handling

-> high precision and robust

Laboratory Automation

-> high flexible, scalable & reliable

Reader Technology

-> high sophisticated and easy to use

CyBi[®]-Screenmachine

CyBi[®]-Lumax

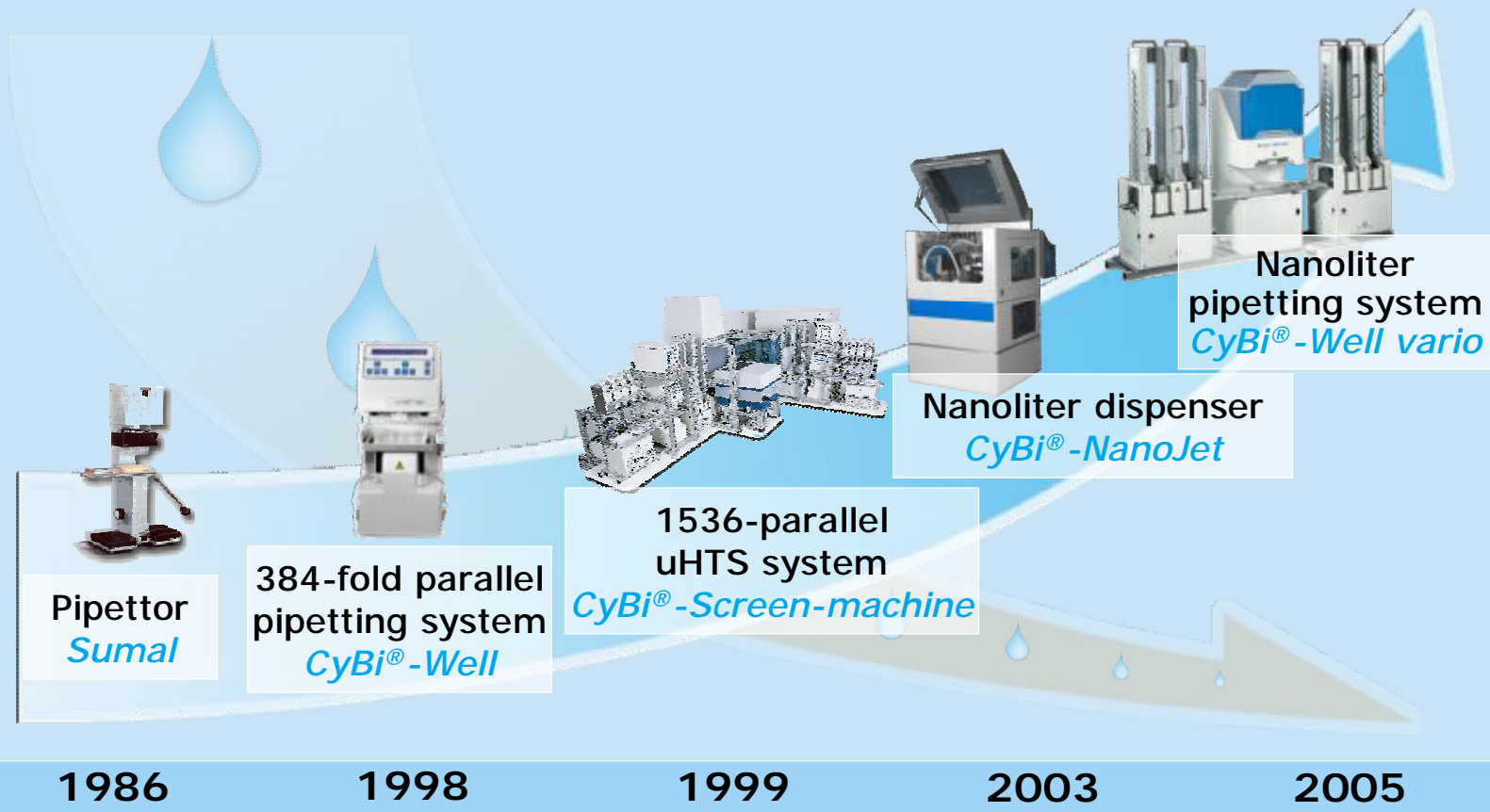
1986


now

Miniaturization in Liquid Handling

more well per plates <- versus -> liquid volume


Solutions for Liquid Handling

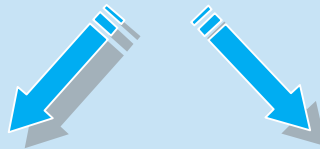Simultaneous pipetting:

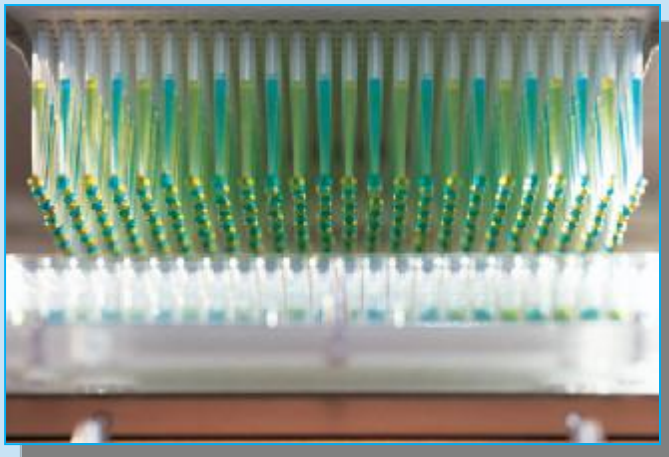CyBi®-DiluS^{pro}
CyBi®-Well
CyBi®-Well vario
CyBi®-Disk

Non-contact dispensing:

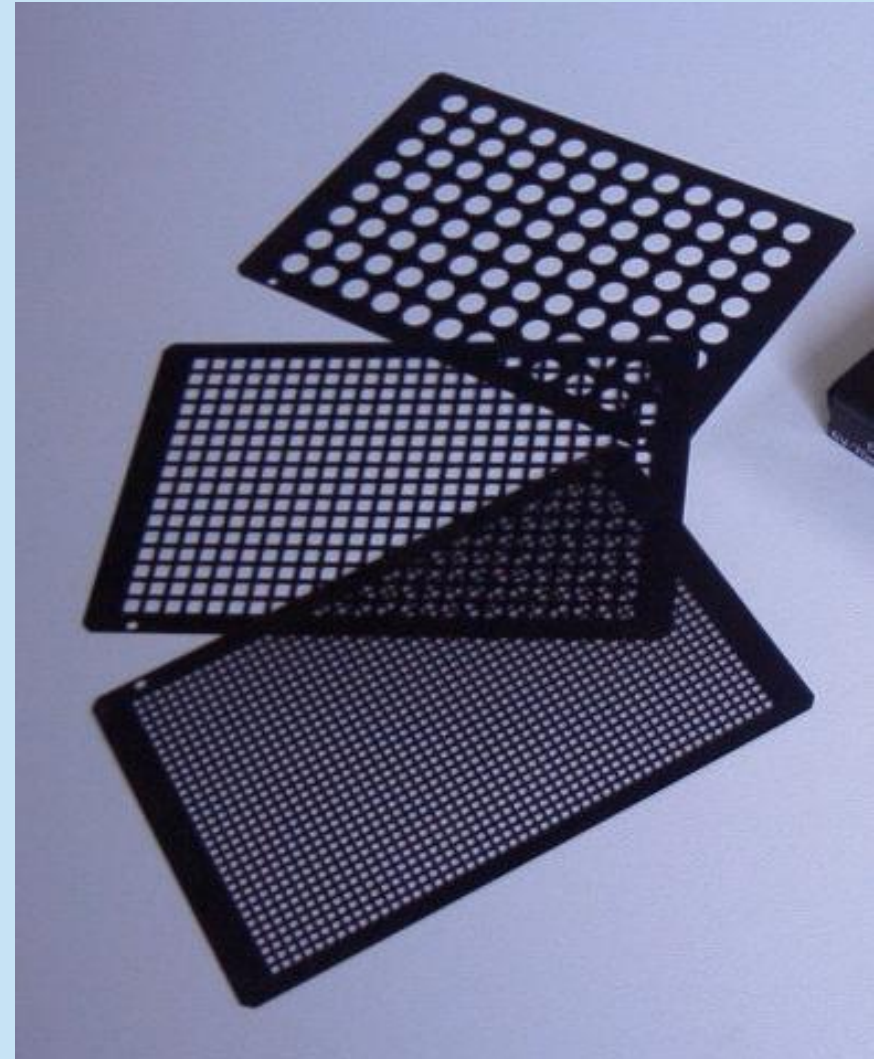
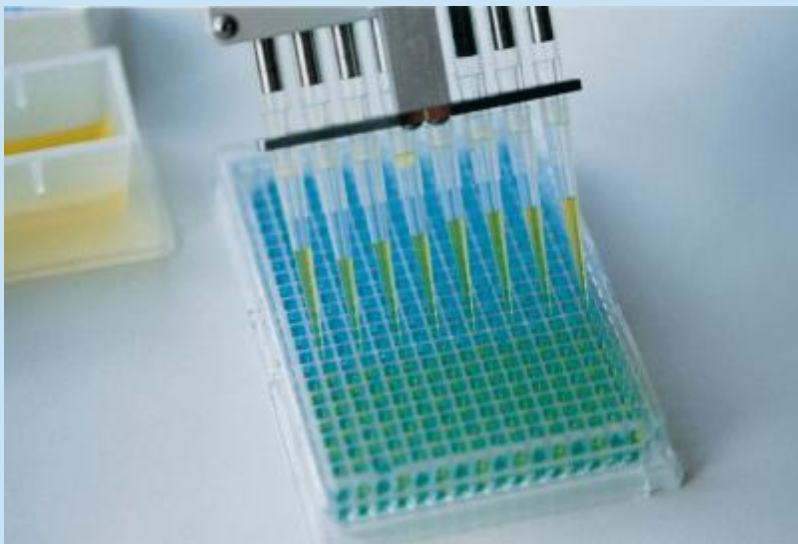
CyBi®-NanoJet
CyBi®-Drop

Liquid Handling - Goals

Fill liquid into wells:



- fast and reliable
- precise und accurate
- find well in 1536 plate
- use disposable cheep tips
- etc...

Liquid Handling - Challenges

Use pipetting or dispensing robotic systems with the goals:

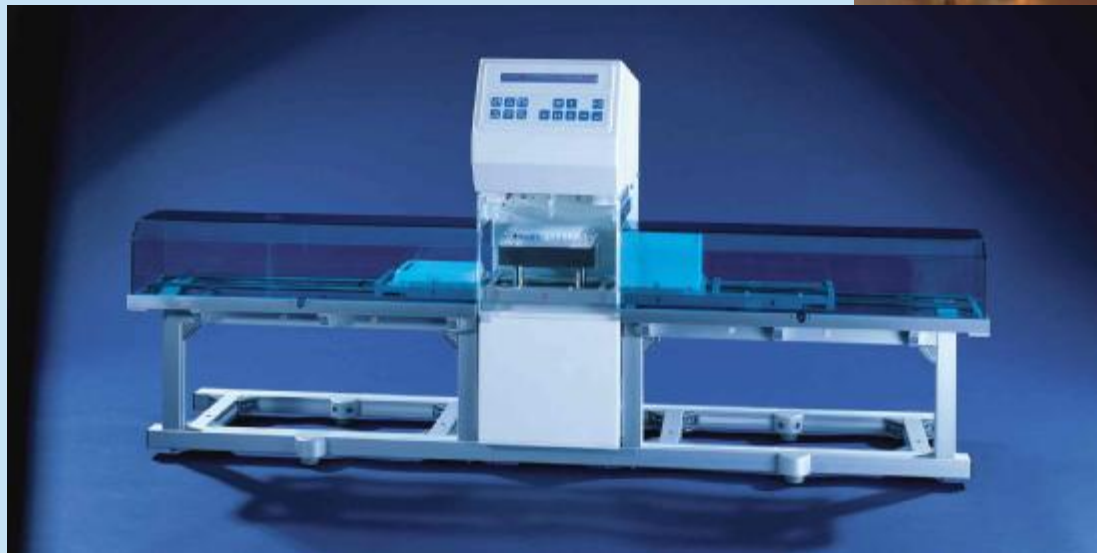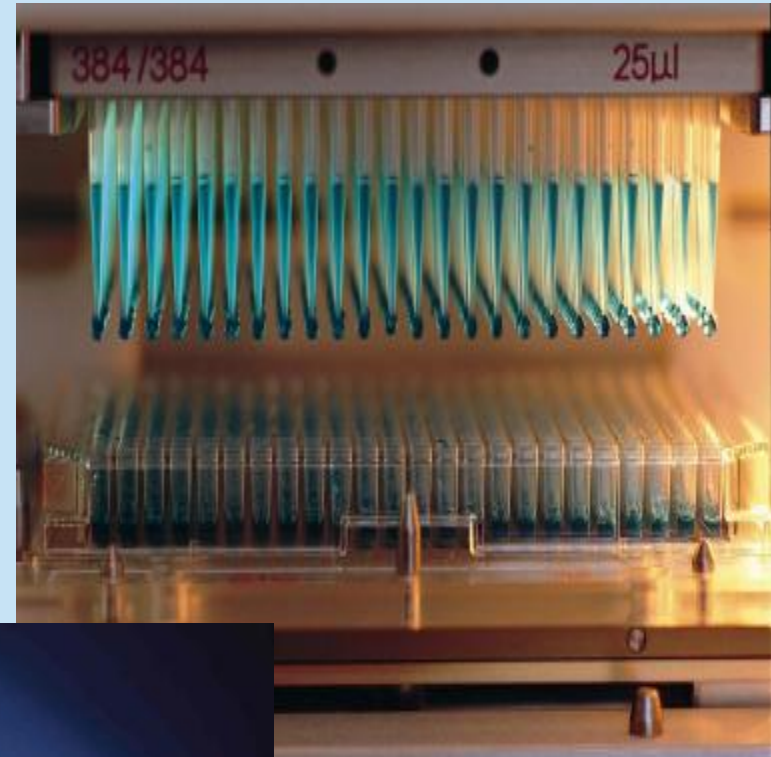- fast and reliable
- precise und accurate
- find well in 1536 plate
- use cheap disposable tips
- etc...



Disposable Tips

Features:

- polypropylene
- injection moulding
- cheep
- one time use

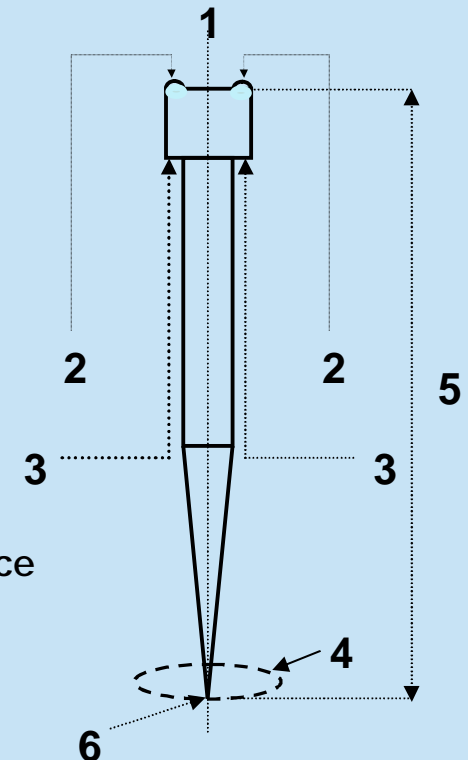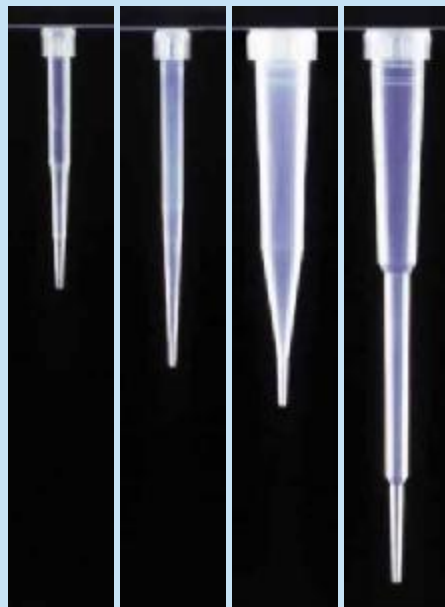- PCR certified
RNAse, DNase and
human DNA free
- sterile
- inline filter


- 1: Straightness
- 2: Evenness of sealing surface
- 3: Evenness of sustaining surface
- 4: Concentricity tolerance
- 5: Length
- 6: No burrs at the outlet


Disposable Tips - Principles

Minimized dead volume between piston and liquid.

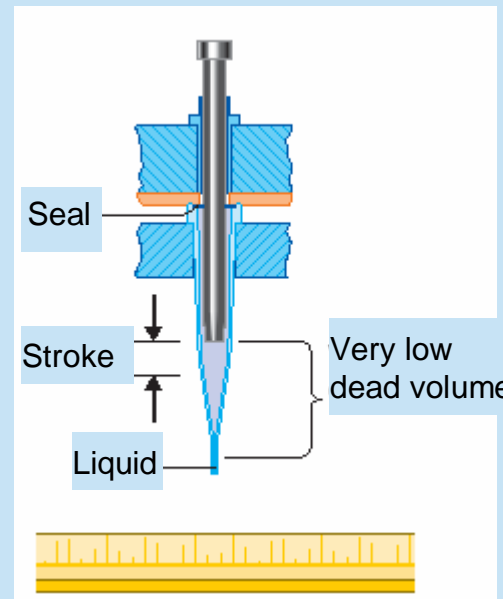- Minimized volume of compressible air combined
- Extremely precise movement of the piston

CyBi®- Well technology

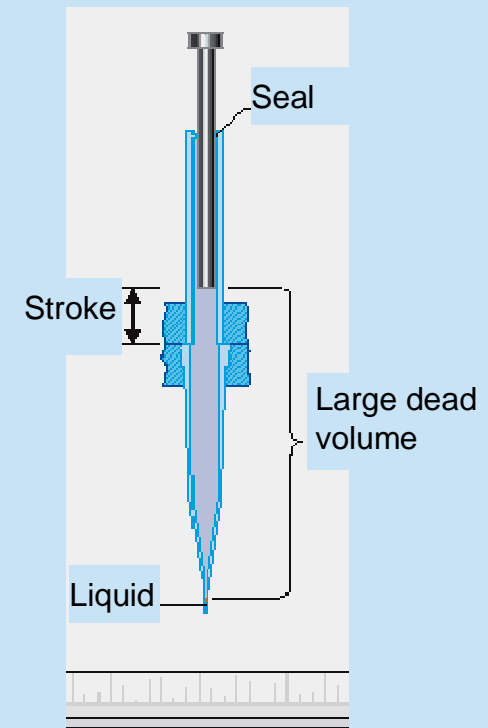Large dead volumes reduce precision

nl-Pipetting – High End Liquid Handling

Nanoliter Pipetting Head: 96/nl head
384/nl head

Volume range: 25 nl – 2500 nl in 1 nl increments

Precision:* 2500 – 500 nl < 3%
500 – 250 nl < 5%
250 – 100 nl < 7%

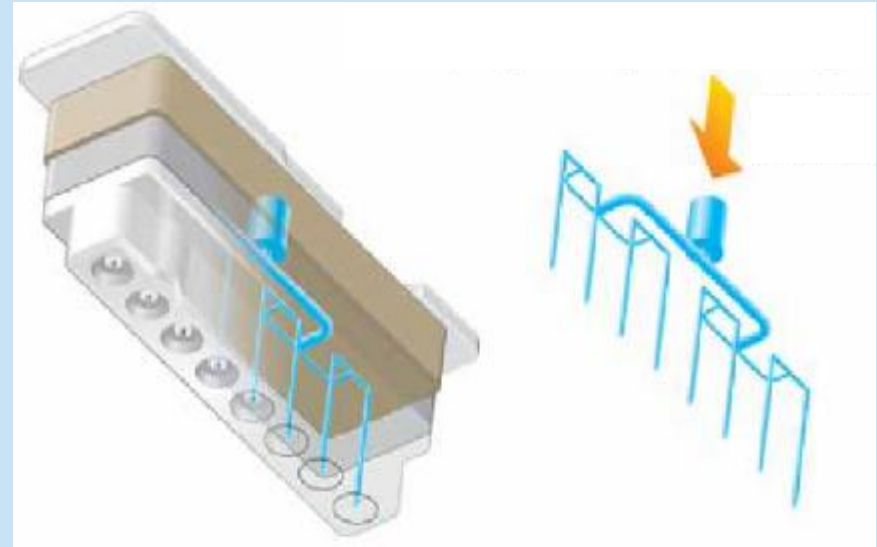
* percentage of standard deviation from the arithmetic mean

Using special ceramic tips

Dispensing Comb

Patented liquid distribution system within the dispensing heads ensures maximum precision and reproducibility



- Dispensing of:
- liquids
 - cell suspension
 - bead suspension

Dispensing Pump

Ceramic rotating piston pump:

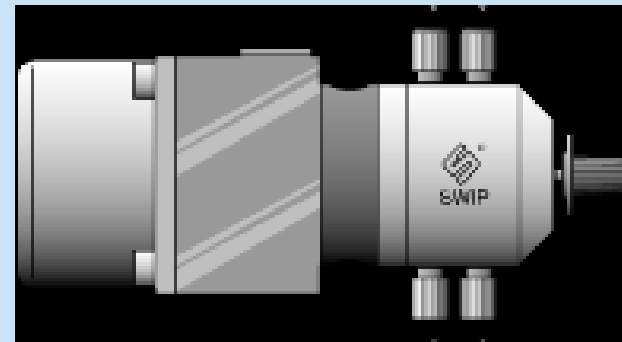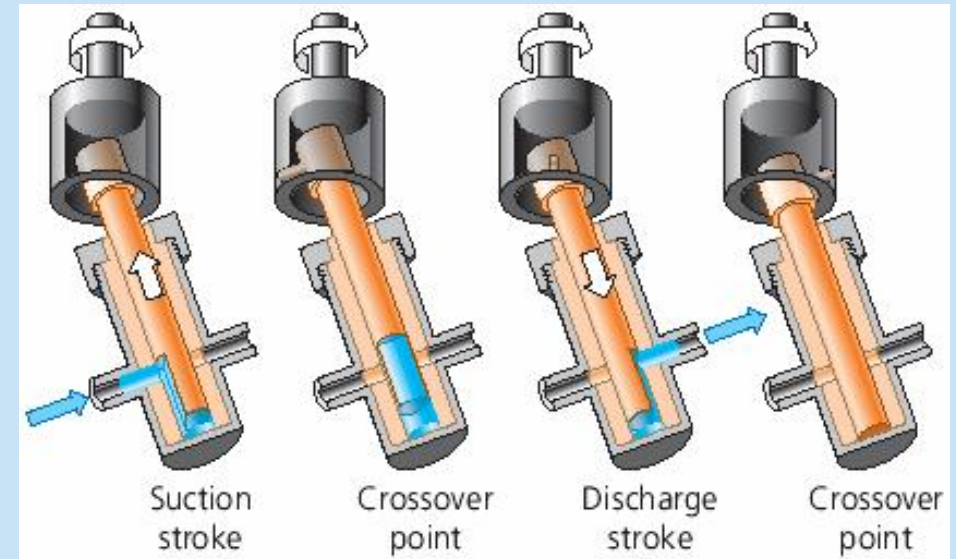- inert
- precise
- robust



Identification

- Identification and tracking of micro plates for
- compound management in automated storage systems
 - interpret assay plates

today : bar code

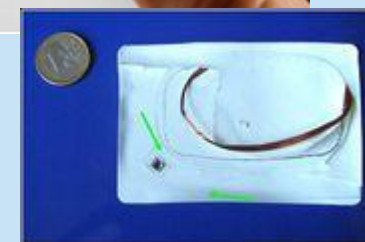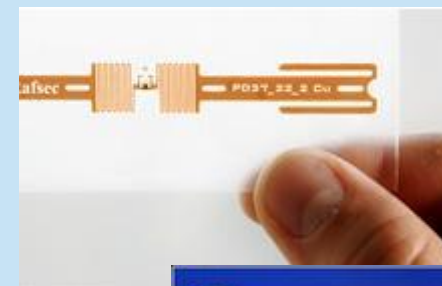future: RFID




storage room with thousands of compounds

Dispensing Technology

Piezo Dispenser

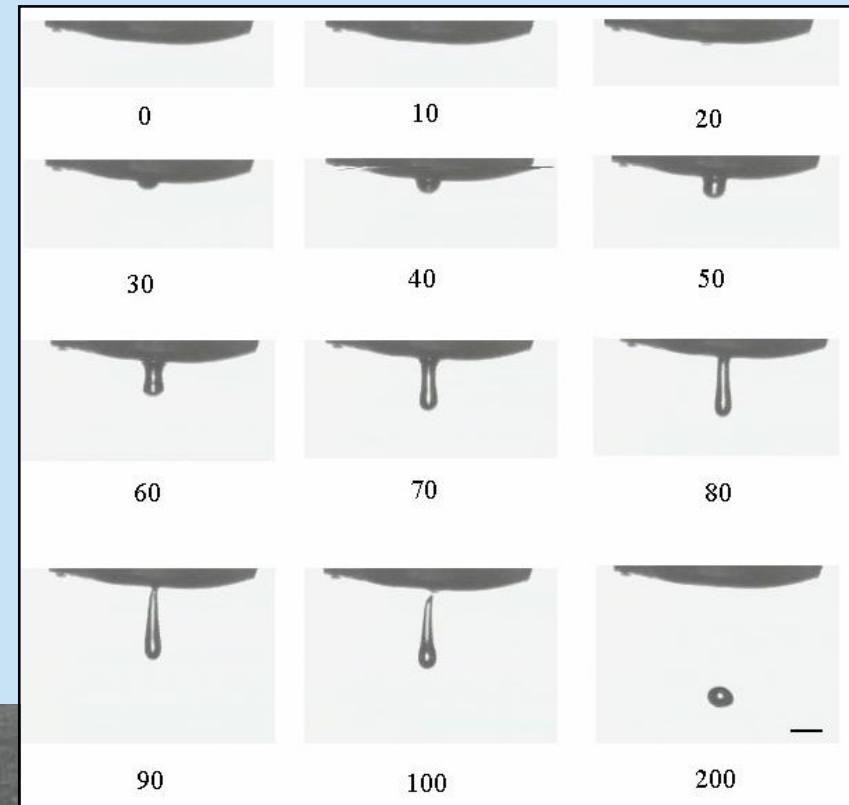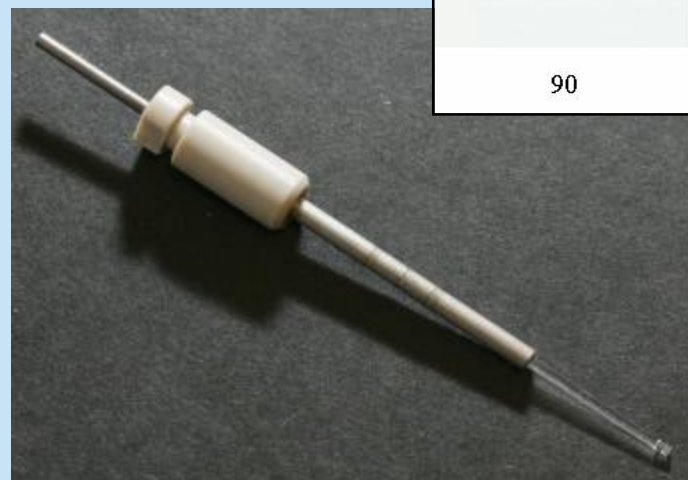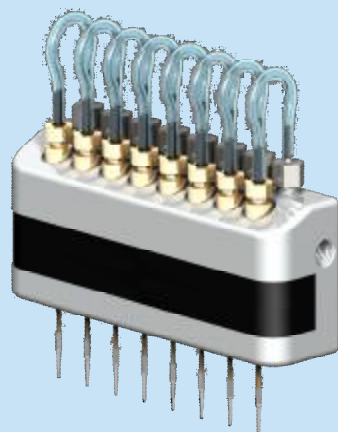- 500pl drops (depends on diameter and piezo force)
- 1nl - 100µl dispensing volume with multi drop
- very fast, very accurate
- needs calibration for every channel
- sometimes unexplainable dropouts

Thermal Measurement

Goal:

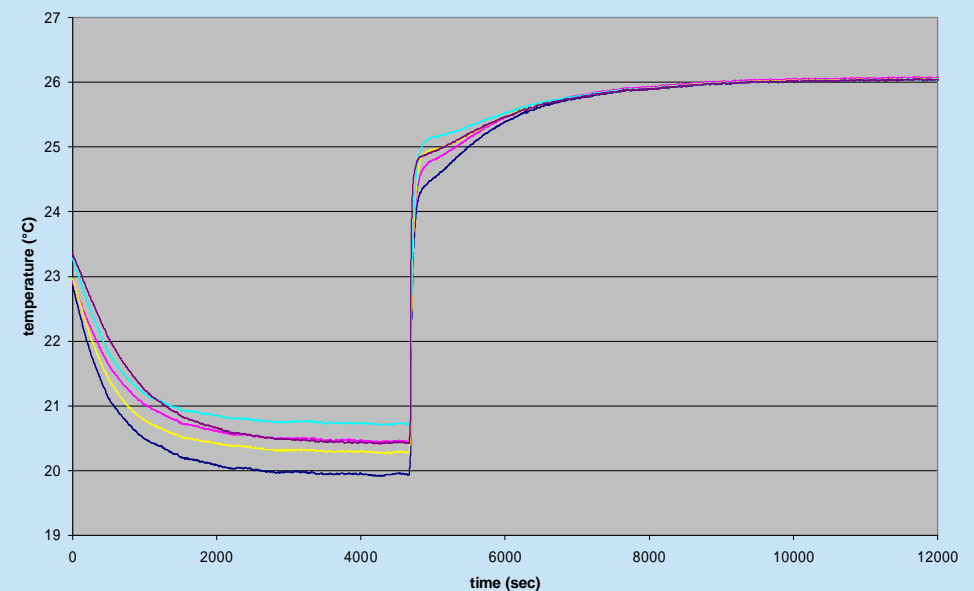
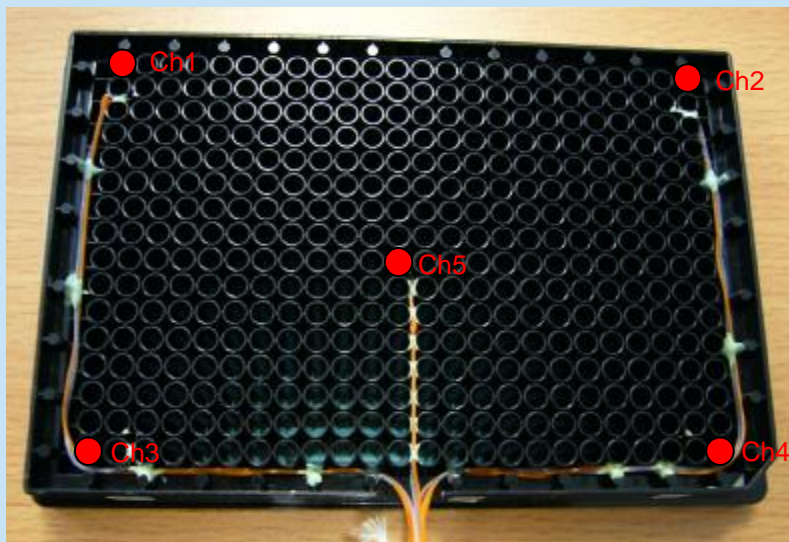
- measurement of gradient in plate
- measurement of equilibration time

Needs:

- fast response
- precise calibrated sensors
- very low thermal capacity
- exactly replaceable

cooling down and equilibration on carousel level 5

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CyBio AG